

Simultaneous Mapping of On- and Off-Target Structural Variants, Insertions and Translocations in Engineered CAR-T Cells

Christopher Tompkins, PhD (Chief Technology Officer, KromaTiD)
 Erin Cross (Senior VP of Research and Quality, KromaTiD)
 Stephen Gross, PhD (Product Research Scientist, KromaTiD)


KromaTiD

Direct, Definitive Genomics

ABSTRACT

Structural variants, such as inversions and translocations, are common by-products of CAR T-cell engineering processes such as CRISPR-Cas9 editing. These variants are formed by misrepair of DNA breaks, for instance between the TRAC loci and an off-target site. These can result in a heterogeneous mixture of low-prevalence variants that involve edit-site, off-target and random breaks. Directional Genomic Hybridization™ (dGH™) is a unique cytogenetic technique for mapping the structure and structural variation of many individual genomes in single cells. Based on images of fluorescently labeled DNA probes designed against a reference genome for specific loci and hybridized to metaphase chromosomes, dGH in-Site™ assays provide true *de novo*, unbiased detection of structural variants or CAR insertions as small as 2kb anywhere in the genome. dGH in-Site™ is ideal for measuring these by-products of editing, mapping the on- and off-target locations of transgenes as well as potentially genotoxic outcomes such as sub-clonal outgrowth, insertion at potentially oncogenic sites and chromothripsis.

directional Genomic Hybridization (dGH™)

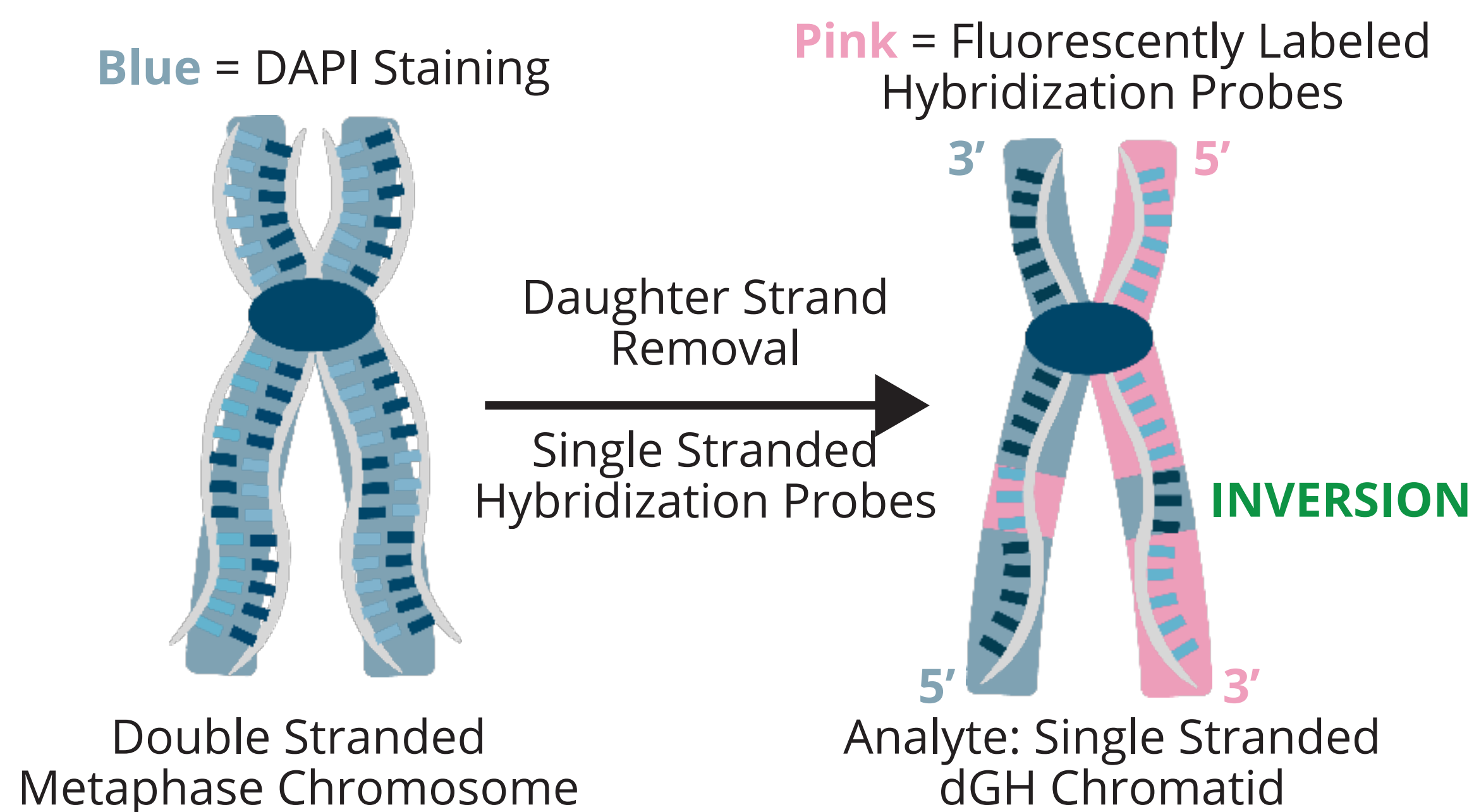

- dGH™ chromosomes contain only the 2 parental DNA strands in opposite orientation.
- Single stranded probes are designed to target only the parental strand (no reverse complements).
- Signals appear on only one chromatid. Inversions present as signal from the opposite sister chromatid.


Comprehensive Analysis Package

	Sequence Variation	Mis-repairs and DNA Damage	Karyotyping
Sequence Based Techniques	NGS (Guide-Seq., Circle-Seq.) ddPCR (Biorad)		
Hybridization Mapping	Genomic Barcoding (Nanostring, Bionano, Genomic Vision)	BAC FISH, SKY (Metaphase, Interphase)	
		PinPoint FISH™ (PPF) (Metaphase, Interphase)	
		dGH in-Site™ (Metaphase)	
		dGH SCREEN™, dGH DSCVR™ (Metaphase)	
Chromosome Staining		G-Banding (Metaphase)	

dGH in-Site™ Strategy

- Sequence:** identify target chromosome location & copy number variation (CNV) by probe color.
 - Loci Marker Probe (Green)
 - Non-genomic Insert (Yellow)
 - Centromere Enumeration Probe (Pink)
 - Telomere Enumeration Probe (Orange)
- Orientation:** opposite sister chromatid indicative of inverted target.
 - TEP signals indicate **BREAKPOINT** in telomeric region. Probe signal on the opposite sister chromatid indicates **INVERSION**.
- Location:** map out Loci positions on chromosome arms.
 - CEP identifies chromosome, position on Q-arm **NORMAL**.

dGH™ Workflow

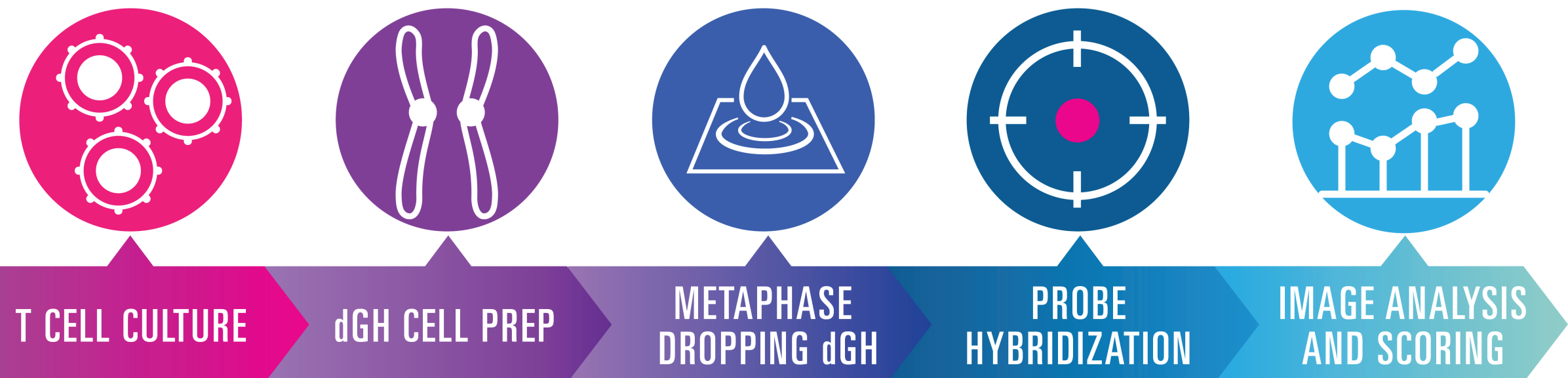

dGH in-Site™ for CAR-T

Verify the Structure of Target Loci + Measure & Locate CAR Insertions = Single-cell Measurements of Many Cells

TRAC: marker probe spans 0.85Mb
B2M: marker probe spans 1.1Mb

- Inversions
- Translocations
- Chromosomal copy number
- Target loci copy number
- On-target insert verification



CAR Transgene Probes

- Insertion signals down to 2kb
- On-target copy number
- Off-target copy number
- Inverted inserts



Percent Occurance of:

- Normal loci
- Inversions/ SCE's of loci
- Translocations of loci
- On-target insertions

Examples: T Cell Metaphase Spreads

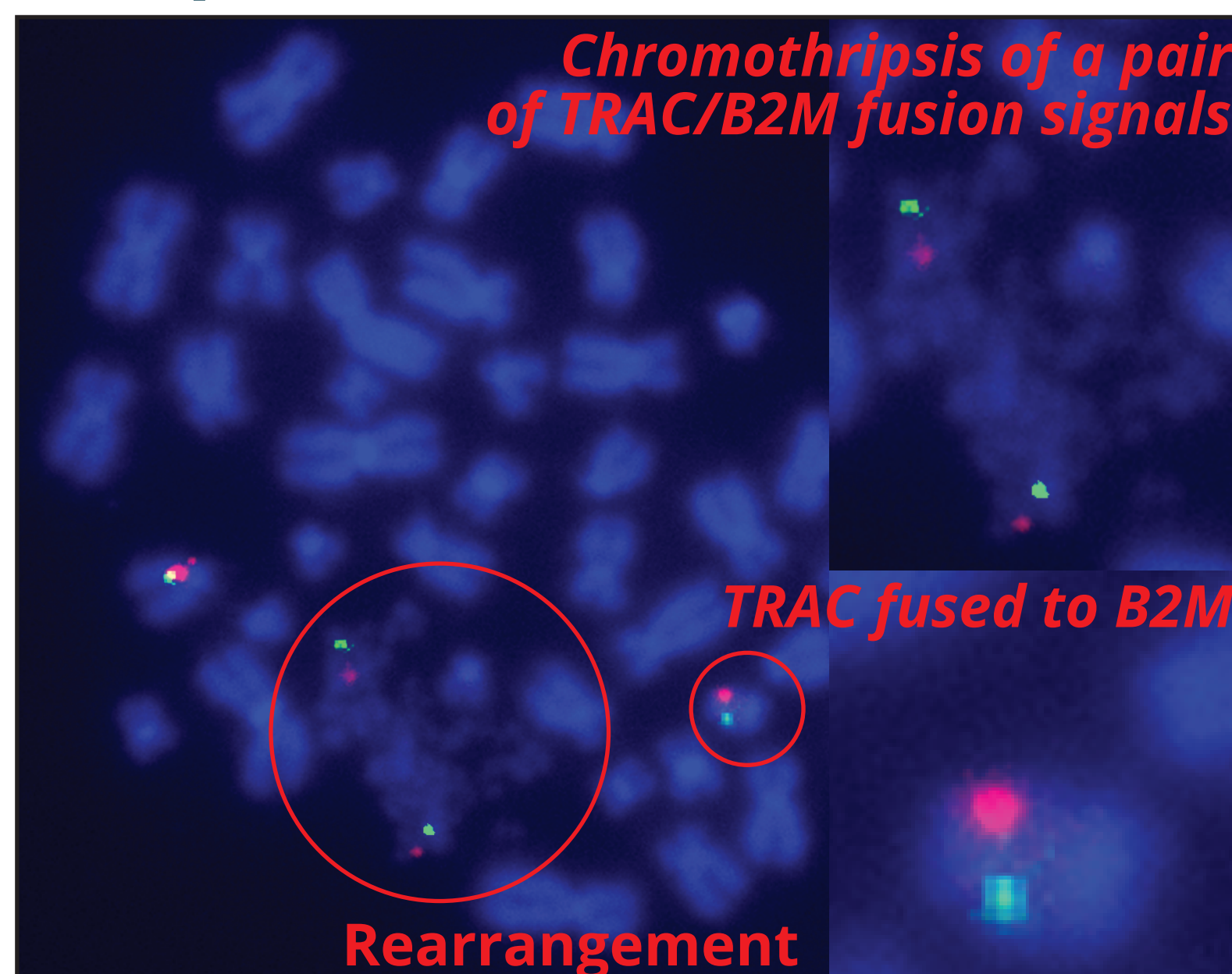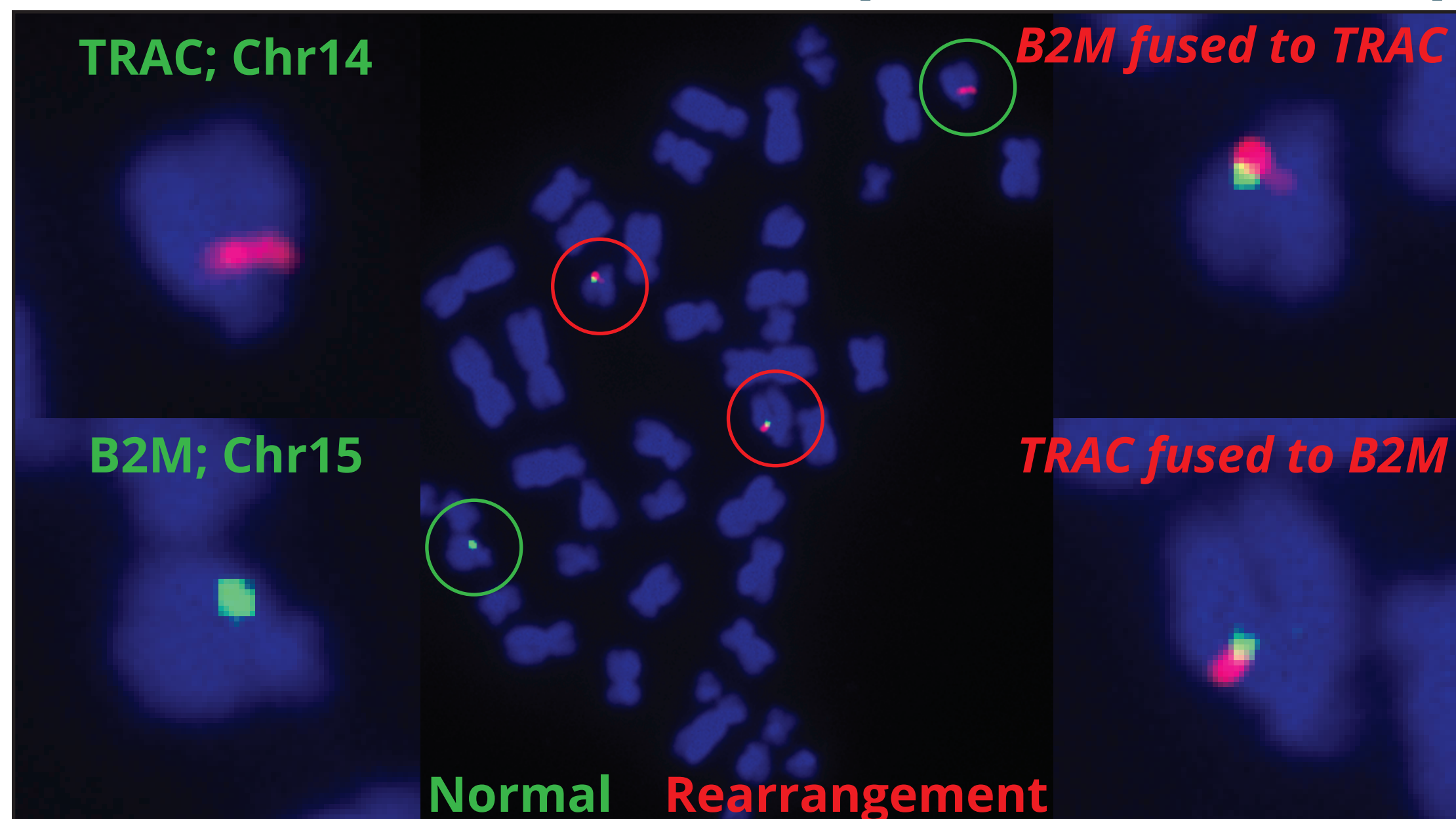

Example: Inserts in iPSCs

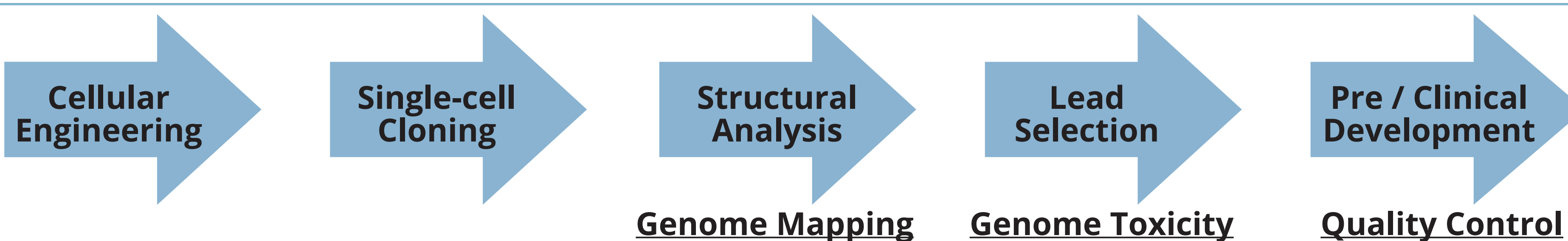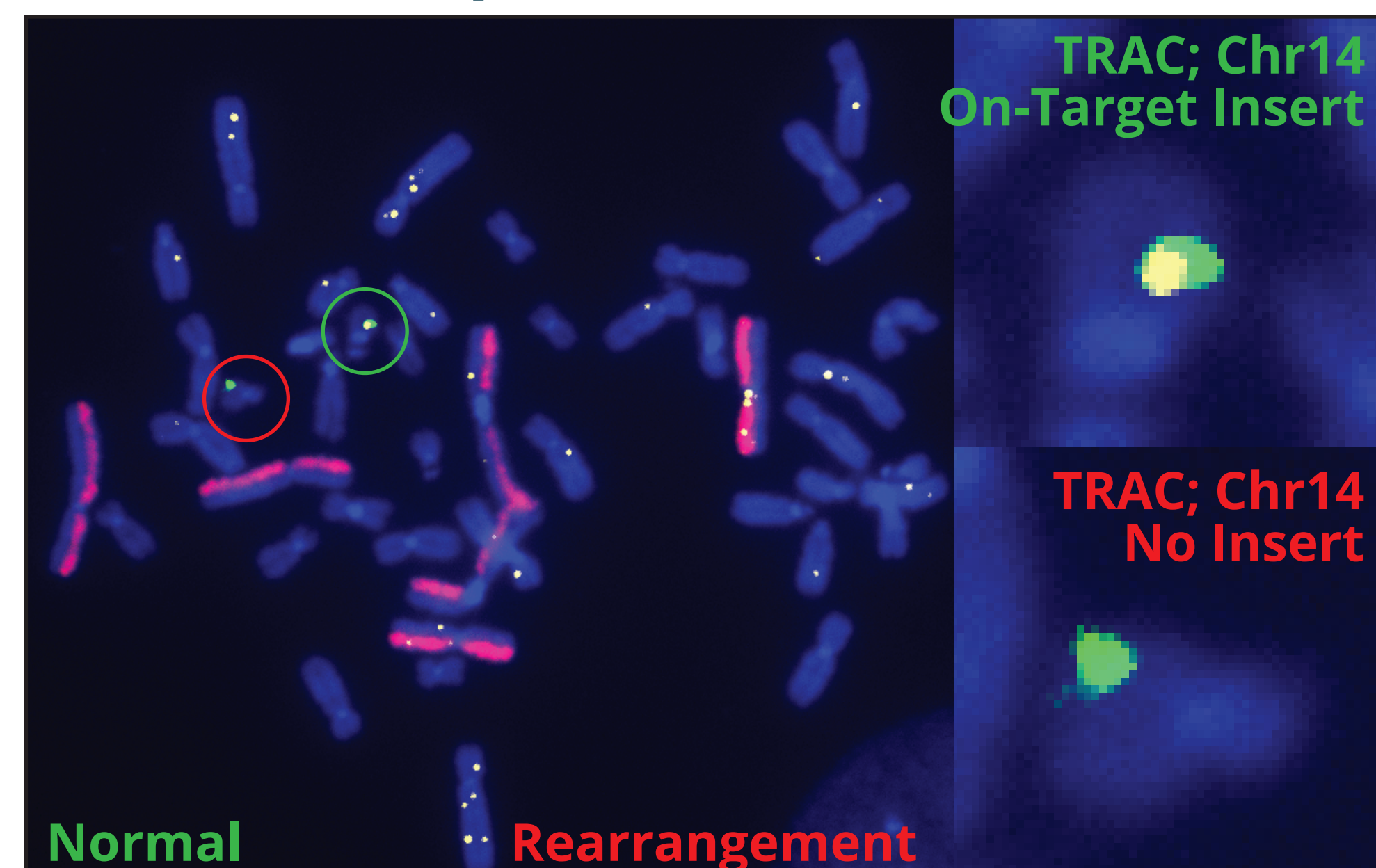

Chromothripsis: Easy to measure with G-Banding, dGH in-Site™ and SCREEN™

Clonal Outgrowth: Measure with dGH in-Site™ timecourse study

Insertional Mutagenesis: Mark high risk loci and track insertions with dGH in-Site™

Structural Variation: Single-cell detection and mapping of variants with dGH in-Site™

Genomic Instability: Early detection of instability with dGH SCREEN™

Aneuploidy: Single-cell detection with dGH SCREEN™ and G-Banding



Please Contact:
 Christopher Tompkins - Chief Technology Officer: ctompkins@kromatid.com
 Dan Vetter - Manager Business Development: dvetter@kromatid.com